

Rules for operating the Innova 4300 shaker:



- 1) This shaker should remain set to 30°C and 300 RPM
- 2) Always use the appropriate sized clamp or rack for the flask or tube that you are shaking. **DO NOT** wedge things in with Styrofoam!
- 3) If there is not an appropriate sized clamp for the flask you want to shake, you can add it. **Be sure to use the right screws for this platform!**
 - a. A tube of the right sized screws can be found taped to the right side of the shaker. The screw should go in easily; if it does not, try a different one.
 - b. Extra clamps and screwdrivers can be found in the cabinet on the south wall of this room.
 - c. If you need a different tube rack, contact Steven Backues and he will put one in for you
- 4) Promptly clean up any spill or splashes or broken glass
- 5) If you are done with the shaker **and** there are no other tubes or flasks in it, you can turn off the shaking (leave the temperature on). But first make sure that there are really no other tubes or flasks in there, or you will ruin someone's experiment.

Contact Steven Backues (sbackues@emich.edu, office 501F) with any questions.